Supel™-Select固相抽出管

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
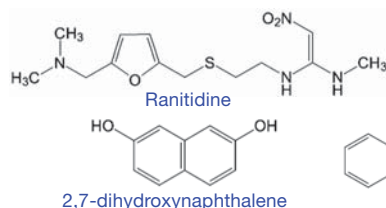
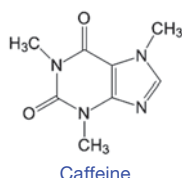
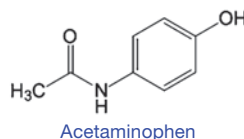
2,7-dihydroxynaphthalene

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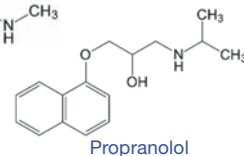Supel-Select HLB

Acetaminophen	98%
Caffeine	96%
Ranitidine	96%
2,7-dihydroxynaphthalene	96%
Propranolol	100%

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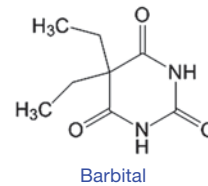
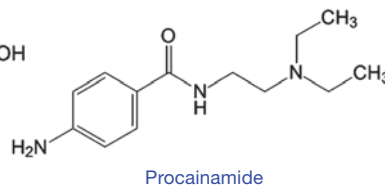
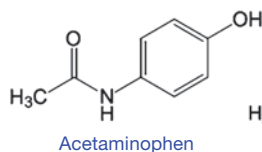
Supel-Select SCX

Acetaminophen	92%
Procainamide	93%
Barbitol	95%

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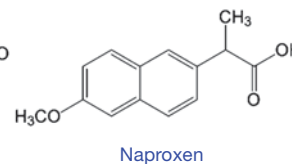
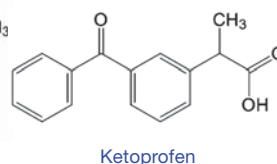
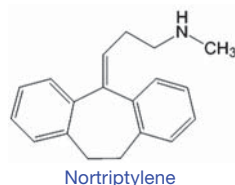
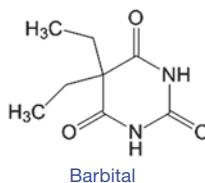
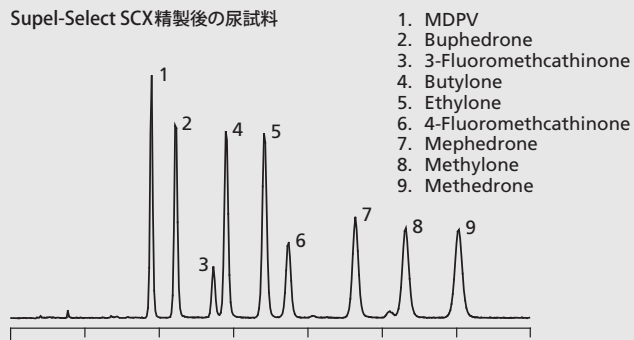
HPLC column : Ascentis® Express HILIC (Si), 10 cm x 2.1 mm I.D., 2.7 µm


Cat.No. : 53939-U

sample/matrix: 1 mL urine spiked to 100 ng/mL of bath salt mixture
 SPE tube: Supel-Select SCX, 30 mg/1 mL (54240-U)
 conditioning: 1 mL 1 % formic acid acetonitrile, then 1 mL water
 sample addition: 1 mL spiked urine
 washing: 1 mL water, 1 mL 1 % formic acid in acetonitrile, 1 mL water
 elution: 2 mL 10 % ammonium hydroxide in acetonitrile
 eluate post-treatment: thoroughly mix via vortex agitation, evaporate 1 mL aliquot to dryness, reconstitute in 500 µL acetonitrile
 column: Ascentis® Express HILIC (Si), 10 cm x 2.1 mm, 2.7 µm (53939-U)
 mobile phase: (A) 5 mM ammonium formate acetonitrile;
 (B) 5 mM ammonium formate water; (98:2, A:B)
 flow rate: 0.6 mL/min
 pressure: 127 bar
 column temp: 35 °C
 detector: MS, ESI(+), 100-1000 m/z
 injection: 1 µL
 sample: 200 ng/mL in acetonitrile

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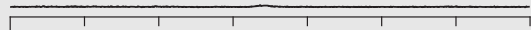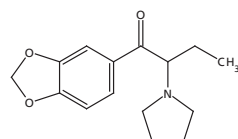
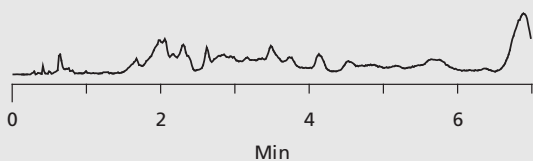
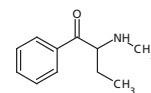
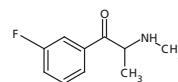
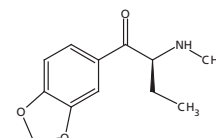
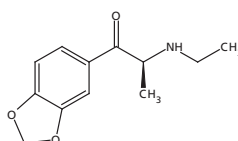
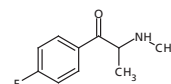
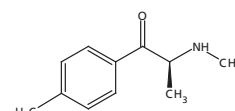
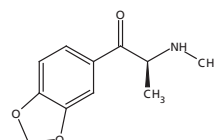
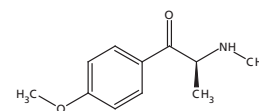
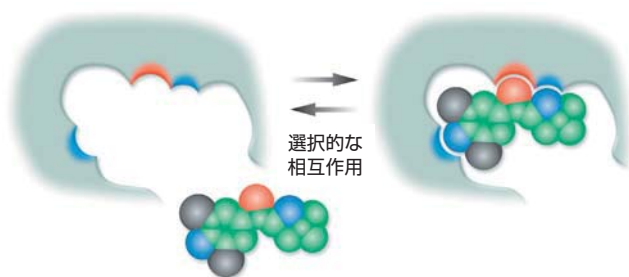
分子鑄型ポリマー (Molecularly Imprinted Polymer)

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© 2014 Sigma-Aldrich Co. LLC. All rights reserved. SIGMA, SAFC, SIGMA-ALDRICH, ALDRICH, and SUPELCO are trademarks of Sigma-Aldrich Co. LLC, registered in the US and other countries. FLUKA is a trademark of Sigma-Aldrich GmbH, registered in the US and other countries. SAFC brand products are sold through Sigma-Aldrich, Inc. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see product information on the Sigma-Aldrich website at sigma-aldrich.com. Ascentis and SupelMIP are registered trademarks of Sigma-Aldrich and Sigma-Aldrich Biotechnology, LP. Supel is a trademarks of Sigma-Aldrich and Sigma-Aldrich Biotechnology, LP.

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